

Laser Power Measurement Spheres



Collecting the
total radiant power
emitted by highly divergent
or collimated sources

Accurate

Independent of beam geometry and without the need for auxiliary optics, Labsphere's Laser Power Measurement Spheres are ideal for collecting the total radiant power emitted by highly divergent sources, such as laser diodes as well as a collimated laser beam. Accurate, repeatable measurements may be performed quickly and easily without painstaking positioning and alignment procedures. Integrating sphere measurements of laser power also eliminate polarization effects often associated with filter attenuation techniques.

Labsphere's Laser Power Measurement Spheres are offered for applications in the wavelength regions from UV-VIS-NIR to the NIR-MIR.

Easy to use

Each laser power measurement sphere is designed to accommodate any standard Labsphere detector assemblies.

The spheres include two statically placed detector ports. Users can mount two detectors simultaneously to measure laser sources with difference in output wavelength. The sphere can also be used for laser characterization by mounting a photodetector on one port and a spectrometer on the other.

A laser target fits into a port opposite the input beam and is designed to receive the first strike from the laser and disperse its energy into the sphere.

Normal tap water or a refrigerated coolant such as ethylene glycol may be used as a coolant. If a recirculator is used, consult the manufacturer as to the choice of coolant. The measurement device is ideal for output measurements from CO₂, Nd:YAG, and other infrared lasers with power density less than 2 kW/cm². The maximum coolant pressure is 15 psi. and the relief valves lifts at 35 psi. Pressure limitation and the coolant outlet temperature should be considered when selecting a coolant base.

Value

- Results independent of beam geometry up to 40° half angle
- Efficiently collects total power from highly divergent sources
- Ability to mount two different detectors or spectrometers simultaneously
- Attenuates the signal to avoid detector saturation

Applications

- Laser and laser diode output characterization
- Optical bench detection

Ordering Information


Model Number	Order Number
3P-LPM-020-IG	AS-02490-001
3P-LPM-020-SF	AS-02489-001
3P-LPM-020-SL	AS-02488-001
3P-LPM-040-IG	AS-02496-001
3P-LPM-040-SF	AS-02492-001
3P-LPM-040-SL	AS-02491-001
3P-LPM-060-IG	AS-02497-001
3P-LPM-060-SF	AS-02495-001
3P-LPM-060-SL	AS-02493-001

Specifications

	LPM-XXX-SL
Sphere Diameters, respectively:	2, 4, 6 inch
Input Aperture Diameters, respectively:	0.5, 1, 1 inch
Sphere Material/Coating:	Spectralon
Optimum Spectral Range:	250 - 2500 nm

LPM-XXX-SL


The LPM-XXX-SL Laser Power Measurement Spheres are integrating spheres designed for the measurement of lasers and laser diodes in the UV-VIS-NIR spectral region (300 - 1800 nm). The interior of the sphere is machined from Spectralon®, a thermoplastic resin with reflectance of approximately 99% over the spectral range of 400 to 1500 nm, and > 95% from 250 to 2500 nm.

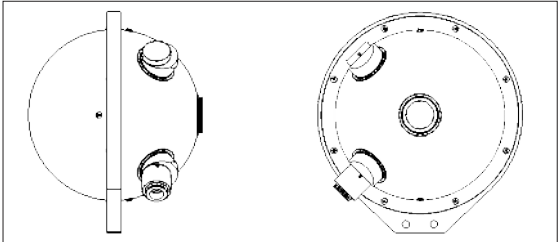

Example of 6 inch LPM

LPM-XXX-SF


The LPM-XXX-SF Laser Power Measurement Spheres are integrating spheres designed for the measurement of lasers and laser diodes in the UV-VIS-NIR spectral region. The interior of the sphere is coated with Spectrafect®, a diffuse white coating applicable over the spectral range of 350 to 1800 nm.

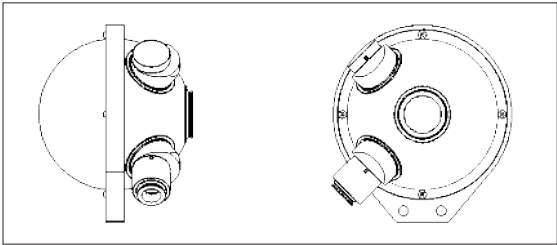

Example of 4 inch LPM

LPM-XXX-IG

The LPM-XXX-IG Laser Power Measurement Sphere is coated with Infragold® diffuse reflective coating, which exhibits high reflectance values (93%–96%) over the NIR-MIR wavelength region.

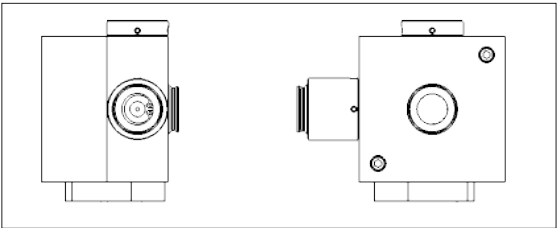

Example of 2 inch LPM

	LPM-XXX-SF	LPM-XXX-IG
Sphere Diameters, respectively:	2, 4, 6 inch	2, 4, 6 inch
Input Aperture Diameters, respectively:	0.5, 1, 1 inch	0.5, 1, 1 inch
Sphere Material/Coating:	Spectrafect	Infragold
Optimum Spectral Range:	350 to 1800 nm	700 nm - 20 μm